

Introduction to the Microscope Lab Activity

Introduction

"Micro" refers to **tiny**, "scope" refers to **view or look** at. Microscopes are tools used to enlarge images of small objects so as they can be studied. The compound light microscope is an instrument containing **two lenses**, which magnifies, and a variety of **knobs to resolve (focus)** the picture. Because it uses more than one lens, it is sometimes called the compound microscope in addition to being referred to as being a light microscope. In this lab, we will learn about the proper use and handling of the microscope.

Instructional Objectives

- Demonstrate the proper procedures used in correctly using the compound light microscope.
- Prepare and use a wet mount.
- Determine the total magnification of the microscope.
- Explain how to properly handle the microscope.
- Describe changes in the field of view and available light when going from low to high power using the compound light microscope.
- Explain why objects must be centered in the field of view before going from low to high power using the compound light microscope.
- Explain how to increase the amount of light when going from low to high power using the compound light microscope.
- Explain the proper procedure for focusing under low and high power using the compound light microscope.

Materials

- Compound microscope
- Glass slides
- Cover slips
- Eye dropper
- Beaker of water
- The letter "e" cut from newsprint
- Scissors

Procedures

I. Microscope Handling

1. **Carry the microscope with both hands** --- one on the arm and the other under the base of the microscope.
2. One person from each group will now go over to the microscope storage area and properly **transport one microscope to your working area.**
3. The other person in the group will **pick up a pair of scissors, newsprint, a slide, and a cover slip.**

Part II. Preparing a wet mount of the letter "e".

1. With your scissors **cut out the letter "e" from the newspaper.**
2. Place it on the **glass slide** so as to look like (e).